

**FACTORY ACCEPTANCE TESTING**  
**DOCUMENTATION**

**PARTICULATE MONITORING SYSTEM**

**XXXXXXXXXXXXXXXXXX**

XXXXXXXXXXXXXXXXXX



**Issue:** Version 1.0

**Date:**

**Reference:** FAT

Developed By	Job Title	Signature	Date

Reviewed By	Job Title	Signature	Date

**Customer Approvals: XXXXXXXXXXXXX**

Approved By:	Job Title:	Signature:	Date:

# Table Of Contents

Item	Description	Page
<b>1</b>	<b>Revision History</b>	<b>4</b>
<b>2</b>	<b>Introduction</b>	<b>4</b>
<b>3</b>	<b>Hardware Configuration Tests</b>	<b>5</b>
3.1	Inspection - Cabinet Build Quality	5
3.2	Inspection - External Layout	6
3.3	Inspection - Internal Layout	7
3.4	Earth Bonding	9
3.5	Signal / Screen Earth Bar	10
3.6	Power Supplies	11
3.6.1	AC Power Supply	11
3.6.2	24 Volt DC Power Supply	12
3.6.3	Redundant Power Supply	14
<b>4</b>	<b>Input / Output Tests</b>	<b>15</b>
4.1	Cabinet Inputs (0mA & 4.00mA)	15
4.2	Cabinet Inputs (12.00mA & 20.0mA)	17
4.3	Accurate Records	19
4.4	Cabinet Outputs (Beacons)	22
4.5	Power Supply Failure Notification	23
4.5.1	Redundant Power Supplies	23
4.5.2	Cabinet Power	23
4.6	Network Communications	24
4.6.1	Cabinet Network Communications	24
4.6.2	Control Node Network Communications	24
4.6.3	Reporting Node Network Communications	25
4.6.4	View Node Network Communications	25
<b>5</b>	<b>Software Configuration Tests</b>	<b>26</b>
5.1	Security Functions	27
5.1.1	OS Log-on (Server)	27
5.1.2	OS Log-on (Control Node)	27
5.1.3	OS Log-on (Reporting Node)	27
5.1.4	OS Log on (View Node)	28
5.1.5	InTouch Log-on	28
5.1.6	Auto Log-off Timeout	29
5.1.7	InTouch Log-on User Levels	29
5.1.8	Change Password	30
5.1.9	Password Length	31
5.1.10	Change Initial Password Assignment	31
5.1.11	Auto Rejection After N Consecutive Password Failures	31
5.1.12	Password Ageing	31
5.1.13	Date & Time Synchronisation to Server	31
5.1.14	Windows OS Vulnerability	31
5.1.15	Audit Log	32
5.1.16	Version Control	32

Item	Description	Page
5.2	Navigation, Operation and Functions - Logged On	33
5.2.1	Control Screen - Logged On	33
5.2.1.1	Suite Overview Screen	36
5.2.1.1.1	LYO Zoom Screen	40
5.2.1.1.2	Aseptic Vial Processing Zoom Screen	43
5.2.1.1.3	Clean Prep and Aseptic Staging Zoom Screen	46
5.2.1.1.4	Formulation & Buffer Prep Zoom Screen	48
5.2.1.1.5	Real Time LYO Screen	51
5.2.1.1.6	Real Time Aseptic Staging & LYO Aisle Screen	53
5.2.1.1.7	Real Time Filling Line & Buffer Table Screen	55
5.2.1.1.8	Real Time Sterilisation Tunnel & Discharge Screen	57
5.2.1.1.9	Real Time Clean Prep Screen	59
5.2.1.1.10	Real Time Formulation & Buffer Prep Screen	61
5.2.1.1.11	Alarm Viewer Screen	63
5.2.1.1.12	Historical Trend LYO Screen	68
5.2.1.1.13	Historical Trend Aseptic Staging & LYO Aisle Screen	78
5.2.1.1.14	Historical Trend Filling Line & Buffer Table Screen	88
5.2.1.1.15	Historical Trend Sterilisation Tunnel & Discharge Screen	98
5.2.1.1.16	Historical Trend Clean Prep Screen	108
5.2.1.1.17	Historical Trend Formulation & Buffer Prep Screen	118
5.2.1.2	System Diagnostic Display Screen	128
5.2.1.3	Windows Controls Screen	129
5.2.1.4	Alarm Point Control Screen	130
5.2.1.5	System Status & Information Panel	133
5.3	Navigation, Operation and Functions (Not Logged On)	150
5.3.1	Control Screen (Not Logged On)	150
5.4	Particle Count Alarms, Displays and Alarm Outputs	152
5.5	Temperature Controlled Counter (Temp Sensor)	171
5.6	Alarm Exception Report	172
5.7	Data Storage	173
5.8	Backup	178
5.9	Software Licences	179
<b>6</b>	<b>Tests Report</b>	<b>180</b>
<b>7</b>	<b>Test Equipment List</b>	<b>182</b>
<b>8</b>	<b>Signatures Record</b>	<b>182</b>
	<u>APPENDIX</u>	
	Copies of Calibration Certificates for Test Instruments	183